



**Procedure**  
**Media Warming Cabinets – Door Catch Repairs**

**Replacement of Door Roller Catch**

Open Cabinet Door

Roll back the door seal in one area and place a flat bladed screwdriver behind the stainless steel door panel

Carefully slide the screwdriver along the door panel, thereby releasing the panel

Completely remove the door panel

Carefully remove the insulation slab material

Unscrew the door roller catch

Fit the new roller catch

Replace the insulation slabs

Slide the door panel into the rubber seal – roll the rubber over the door panel

Check that the door closes correctly onto the stud; adjust stud if necessary (see below)

**Replacement of the Door Stud**

Open the cabinet door

Slacken the stud retaining nut and then unscrew the stud

Screw in the new stud and adjust it so that the door shuts tightly

Tighten the retaining nut

**Please Note**

Warming Cabinet Repairs should only be carried out by suitably qualified Service Engineers